

April 6, 2022

RE: Release of BG Analytics[®] software update Version 1.1.1.57378



BG Analytics[®] is a unique, easy-to-use, intuitive software application that provides a qualitative measurement of (1→3)-β-D-Glucan for use with the Fungitell STAT[®] kinetic assay. BG Analytics[®] software collects and processes data from the Lab Kinetics 8-well incubating tube reader, stores the information in a database and produces reports of the test results.

BG Analytics[®] Version 1.1.1.57378 Release Notes

- Made necessary code changes to address application crashes when installed on a computer with an Intel HID Sensor Collection V2
- Updated the lower cutoff limit for results that were previously below range ending the high frequency of invalid results for low signal Negative clinical samples
- Changed the textual description in the Index field from NaN to “Index Not Calculated”

BG Analytics[®] is now a digital distributed only software application and can be downloaded from the Associates of Cape Cod, Inc., (ACC) software portal on the ACC website at <https://portal.acciusa.com>.

We appreciate your business and will continue to provide you with industry leading products and services to meet your current and future needs.

Sincerely,



Celeste P. Jones
Clinical Diagnostic Product Manager